



## automatic stirrer, magnetic mixer

Model: **APEX-JCX588**

Standard: **GB/T25044**

Introduce: APEX-JCX588 widely used in major universities, environmental protection, scientific research, health and epidemic prevention station, petroleum, chemi...

### Detailed product information

#### APEX-JCX588 Magnetic Mixer

First, automatic stirrer, magnetic mixer use

Magnetic mixer is commonly used [food lab equipment](#). APEX-JCX588 widely used in major universities, environmental protection, scientific research, health and epidemic prevention station, petroleum, chemical, medical and other units, is an ideal laboratory personnel necessary for modern laboratory instruments.

Second, the automatic stirrer, magnetic mixer characteristics

Magnetic mixer is commonly used [food lab equipment](#). APEX-JCX588 magnetic stirrer made of high quality DC motor, low noise, smooth speed control, fully enclosed heating plate can be used for auxiliary heating. Temperature magnetic stirrer can set the temperature and temperature display, long-term heating to use, intuitive and accurate digital display. PTFE and high-quality refined into a stirring magnet, high temperature, wear resistance, chemical corrosion, strong magnetism. Can be adjusted in a closed mixing container, easy to use and very good.

Third, the automatic stirrer, magnetic mixer technical parameters

Power supply: AC220V  $\pm$  10

Stirring speed: 100 ~ 2000r/min

Maximum mixing capacity: 2000ml

Dimensions: 120  $\times$  120  $\times$  65 (mm)

Motor power: 25w, heating power: 150w

Speed: 0 ~ 2400 rev / min, infinitely adjustable.

78-1 magnetic stirrer Description:

Fourth, use of methods

Magnetic mixer is commonly used [food lab equipment](#). When using the APEX-JCX588 magnetic heating stirrer, first check the accessories are complete, then the order first installed fixtures, to the stirred solution needed to put chrome plate in the middle. Added to the solution in the beaker and stir the solution, and then plug in the instrument first plug the power cord, and then turn the power speed power switch, indicator light, began to work, and then select your desired temperature, speed gradually raised by the low-speed, high-speed file directly allowed to start in order to avoid stirring of sync, causing the beating, not heated without stirring, heating has the band switch control, cut off the power when not working. The instrument should be kept clean and dry, non-solution into the machine.

Fifth, automatic stirrer, magnetic mixer notes

(1) found stir stir stir beating or not, check the beaker is stable, the location is positive.

(2) medium speed for extended life

Please connect the ground wire (3) the use of

Sixth, the whole list

1, machine 1 sets

2, a stirring

3, power cord 1

4, a chrome pole

5, chrome ten bytes of a

6, chuck a rubber

7, bakelite support screws 2